The packed column for HPLC

Develosil XG-C1

It is the column which made C1 (Trimethyl) group modify. There is performance which is not in C1 conventional column, and it contributes to many analysis systems.

1. Specification and performance

Develosil XG-C1 has introduced C1 (Tori methyl radical) into the inhouse high grade silica gel.

(USP column code: Equivalent to L13) The retention and separation which cannot be predicted in C1 conventional column by the height of surface area and special end cap treatment are attained.

It corresponds to many analysis kinds, such as a basic compound, a vitamin, and drugs, including an organic acid.

Characteristic of Develosil XG-C1


Surface area	450 m²/g
Pore diameter	10 nm
Pore volume	1.15 mL/g
Ligand	Trimethyl
Carbon content	7.50%
Endcapping	\bigcirc
pH range	рН 1.5 - 8
Use upper limit pressure	20 Mpa
Use maximum temperature	50 ° C

2. The trait of Develosil XG-CN

_	Product Name	Surface area (m2/q)	Pore diameter(nm)	Pore capacity (mg/ml)	Carbon content (%)	Endcap
	XG-C1-5	450	10	1.15	7.3	Specially
L	TMS-UG-5	300	14	1.05	4.5	Double
	TMS-5	350	12	1.05	4	Single

3. STD separation comparison

Develosil C1 column Comparison of standard chromatogram

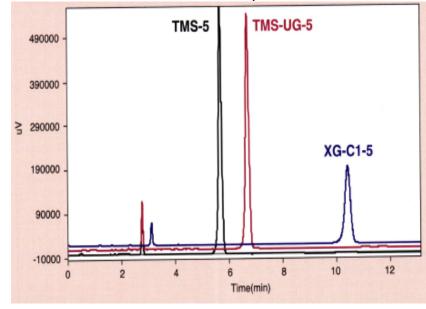


Conditions;		
Column	: Develosil* T	MS-5, TMS-UG-5, XG-C1-5
Size	: 4.6x150mm	
Mobile phase	: Acetonitrile/	Water=70/30
Flow rate	: 1.0ml/min	
Temperature	: 30°C	
Detection	: UV254nm	
Sample	: 1.Benzene	2.1,3,5-Triphenylbenzene

Separation comparison by a STD sample Though it has the same ligand, separation patterns differ, respectively.

If it is XG-C1, about 1.5 times as many retention is shown rather than C1 conventional column. Compared with C1 conventional column, a high grade silica gel with large surface area is adopted as a base material.

Since surface area is large, retention is a column which has the feature of not only a big tree but XG series.

4. Separation characteristics by C1 modification type Packing material

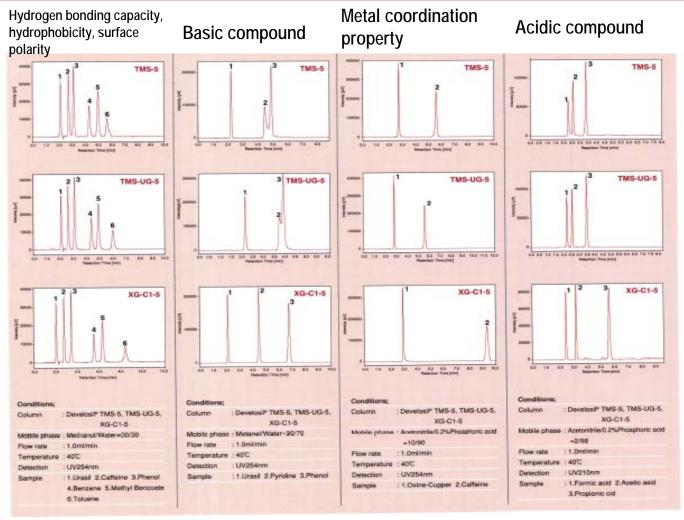


Develosil XG-C1 shows good separation as compared with C1 conventional column.

In separation of a "basic compound", what was insufficiency of separation is separated finely, and the peak shape of pyridine is also sharp. And, the very good result is shown also in separation of an "acidic compound", and the characteristics which are not in C18 or C30 are shown.

X It is carrying out on specific terms. It becomes separable by changing terms also in C1 conventional column.


5. Application

1. Analysis of a pyridine derivative

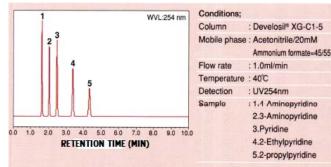

Form also with sharp pyridine represented by the basic compound

3. Analysis of ascorbic acid in a drink, and citric acid

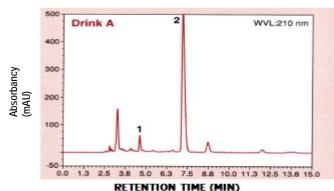


Suppose Sup

Ascorbic acid and citric acid contained in two sorts of commercial item drinks were analyzed.


Inj

4. Analysis of an aromatic compound 5. Analysis of water soluble vitamin

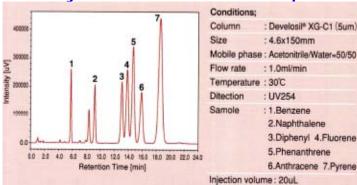


Analysis of various aromatic compounds is possible. All the degrees of separation attain above one.

6. Analysis of drugs

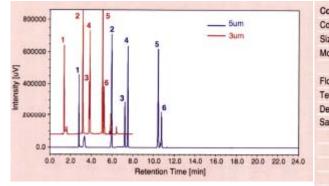


Drugs are analyzed by gradient elution. By using particle diameter 3um, analytical time is shortened sharply.

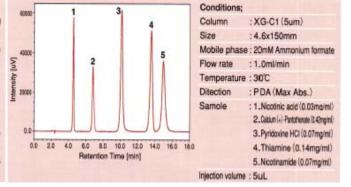

The vitamin analyzed until now using C18 or C30 can be analyzed also by XG-C1.

onditions;				
olumn	: Develosil* XG-C1 (5um, 3um)	Gradient :		
ize	: 4.6x150mm (5um) , 4.6x75mm (3um)	5um		
lobile phase	: A) Water + 0.1%Phosphoric acid	Time (min)	A (%)	B(%)
	B) Acetonitrile + 0.1%Phosphoric acid	0	98	2
low rate	: 1.0ml/min	10	2	98
emperature	: 30°C	15	2	98
etaction	: PDA (Max abs.)	15.1	98	2
ample	: 1.Uracil (0.08mg/ml)			
	2.Acetaminophen (0.10mg/ml)	3um		
	3.Guaifenesin (0.08mg/ml)	Time (min)	A (%)	B(%)
	4.Diphenhydramine (0.10mg/ml)	0	98	2
	5.Flurbiprofen (0.08mg/ml)	3.3	2	98
	6.lbprofen (0.08mg/ml)	4.95	2	98
jection volume	e : 5.0uL	5	98	2

2. Analysis of an organic acid

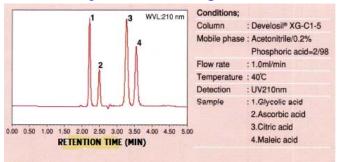

In C18, the organic acid which is hard to separate can also be divided finely.

6. Price List

I.D.2.0mm,3.0mm,4.0mm same price

Convention columns

■ Conventiona columns

Packing name	ing name Column Size		P/N	Price (yen)	
	I.D. 4.6 x 35mm		XGC1346035W	31,000	
	I.D. 4.6 x 50mm		XGC1346050W	33,000	
	I.D. 4.6 x 75mm	ე	XGC1346075W	39,500	
	I.D. 4.6 x 100mm	3	XGC1346100W	46,000	
	I.D. 4.6 x 150mm		XGC1346150W	51,000	
	I.D. 4.6 x 250mm		XGC1346250W	65,000	
XG-C1	I.D. 4.6 x 35mm		XGC1546035W	28,000	
	I.D. 4.6 x 50mm		XGC1546050W	30,000	
	I.D. 4.6 x 75mm	5	XGC1546075W	36,500	
	I.D. 4.6 x 100mm		XGC1546100W	43,000	
	I.D. 4.6 x 150mm		XGC1546150W	48,000	
	I.D. 4.6 x 250mm		XGC1546250W	60,000	

• Guard column for conventiona

Product name	P/N	Price (yen)			
Guard cartridge (4pieces) 4.0x10mm	XGC1540010C	18,000			
Guard cartridge Holder 4.0x10mm	HO00040010C	9,500			
Guard cartridge set 💥 XGC1540010W					
«Guard cartridge set = Holder (1piece) + cartridge 4.0x 10mm (1piece)					

Semi-micro columns

■ Semi-micro columns

Packing name	Column Size	Particle Size	P/N	Price (yen)
	I.D. 2.0 x 35mm		XGC1320035W	31,000
	I.D. 2.0 x 50mm	ŀ	XGC1320050W	33,000
	I.D. 2.0 x 75mm		XGC1320075W	39,500
	I.D. 2.0 x 100mm	3	XGC1320100W	46,000
	I.D. 2.0 x 150mm		XGC1320150W	51,000
	I.D. 2.0 x 250mm		XGC1320250W	65,000
XG-C1	I.D. 2.0 x 35mm		XGC1520035W	28,000
	I.D. 2.0 x 50mm		XGC1520050W	30,000
	I.D. 2.0 x 75mm	5	XGC1520075W	36,500
	I.D. 2.0 x 100mm		XGC1520100W	43,000
	I.D. 2.0 x 150mm		XGC1520150W	48,000
	I.D. 2.0 x 250mm		XGC1520250W	60,000

• Guard column for Semi-micro

Product name	P/N	Price (yen)		
Guard cartridge (4pieces) 1.5x10mm	XGC1515010C	20,000		
Guard cartridge Holder 1.5x10mm	HO00015010C	15,000		
Guard cartridge set X XGC1515010W				
«Guard cartridge set = Holder (1piece) + cartridge 4.0x 10mm (1piece)				

Prep columns

Prep columns

• Prep guard column

Packing name	Column Size	Particle Size	P/N	Price (yen)	Packing name	Column Size	Particle Size	P/N	Price (yen)
XG-C1	I.D. 10.0 x 150mm	5	XGC15P1150W	120,000		I.D. 8.0 x 10mm)mm) 5	XGC1580010W	20,000
	I.D. 10.0 x 250mm		XGC15P1250W	140,000		(for I.D. 10.0mm)			
	I.D. 20.0 x 150mm		XGC15P2150W	245,000		I.D. 20.0 x 50mm		XGC15P2050W	115,000
	I.D. 20.0 x 150mm		XGC15P2250W	300,000		(for I.D. 20.0mm)		X00101 2000W	113,000